


Mounting of Proximity switch on Keofitt valve-head type W9n

1:/ Push the proximity switch into one of the two tracks.



2:/ Turn the little lock into horizontal position and tighten the screw.

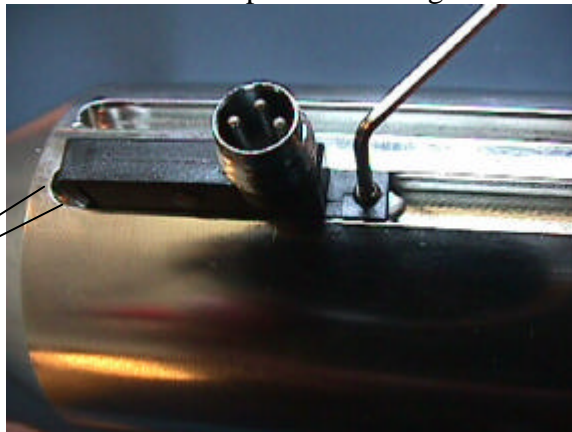

Distance between
prox. switch and
the bottom of the
track app. 5-10 mm

3:/ To test the functionality there is a small light diode on the side of the proximity switch. When the power supply is turn on the diode will light if mounted correctly. By use of the Keofitt lever, open the valve, and the diode will stop lighten. If it doesn't turn off, adjust the proximity switch and try again



Light diode

For electrical connection, see documentation.